

Installing the Spectrum Excel Add-in In Excel 2007

Microsoft Excel 2007 is different from previous versions and requires that you manually add the Spectrum Excel Add-in after installing the software.

Follow these steps and you will be able to access the Spectrum Menu in Excel 2007:

1. Install the Spectrum Excel Add-in software
2. Open Excel 2007
3. Click on the MS Office button located in the top left corner of the Excel window.
A new window opens
4. Click “Excel Options” near the bottom right corner of the window. A new window opens
5. Click “Add-ins” in the left panel of the window
6. Choose “Excel Add-ins” from the drop-list at the bottom of the window and click Go. A new window opens
7. Click **Browse** and browse to C:\Program Files\Veriteq Instruments\SpectrumExcelAddin. Double-click the SpectrumExcelAddin.xla file that is in this directory
8. “Veriteq Spectrum Excel Addin v1.10” now appears in the Add-ins window. Confirm that there is a checkmark beside it click **OK**
9. The Spectrum menu has been added to the Add-Ins tab in the Excel workbook